

REI SYSTEM (Rotor Easy to Install)

Our new REI (Rotor Easy to Install) system for quick and easy exchange of rotors, allows the rotor to be installed and locked securely without the need for tools, as well as unlocked by simply removing it from its position.

Thanks to its ergonomic design, it can be operated comfortably with one hand. The rotor can be removed and replaced in just a few seconds, without rotating the hand, or pressing buttons. It just needs to be positioned on the motor axis and it will be automatically anchored thanks to the REI system.

For greater security, the equipment screen warns the user if the rotor is not properly installed.

To remove it, gently lift the red handle, which will release the REI system allowing the rotor to be removed.

